

CAP TOOL GUIDE

Cap Tool Quick Use Guide

The all new BRAND Cap Tool simplifies the process for attaching and detaching caps on PCR plates. Made of durable polyamide (PA) material, this handy tool has an end specifically designed to seal cap strips onto BRAND standard profile PCR plates, as well as PCR strip caps and plates from other manufacturers. The opposite end features a pointed hook for efficient removal of PCR caps.



Cap sealing end

Cap removal end



Instructions for Use

1. Fill standard profile BRAND PCR plate with experimental samples and reagents.
2. **For PCR cap sealing:** Place PCR caps onto the wells. Using the sealing end, place the Cap Tool directly on top of the cap and apply a slight downward pressure to secure the cap in the PCR well. Repeat the process for all wells containing PCR samples.
3. **For PCR cap removal:** To remove the PCR cap strip from the PCR plate, place the pointed end underneath the strip in between two caps and apply a slight upward pressure. The cap should be released from the PCR plate. Repeat for each well with a cap.



Lab Rats

Trust BrandTech®!

BRANDTECH®
SCIENTIFIC, INC.

Toll Free: 888-522-2726
www.brandtech.com